

Automet[™] Service Plans



Protect Your Investment

Labs run more efficiently and consistently with well-maintained equipment.



Reduce Unplanned Downtime

Regular equipment maintenance will reduce the hassle of unexpected downtime and associated costs.



Ensure Consistent Results

Confidently analyze lab results with the equipment maintained by the trusted experts at Buehler.





Service Plan Options

Identify the best Service Plan based on the lab's needs.

	Preventative Maintenance	Extended Warranty
Visits Per Year	1	2
Wear Parts Kit	\checkmark	\checkmark
Unlimited Phone Support	\checkmark	\checkmark
Replacement Part Discount	10%	25%
Unlimited Labor Repairs		\checkmark
Priority Response		√

- Preventative Maintenance ensures consistent results
- Visit includes Wear Parts Kit and 12 Point and 14 Point Maintenance Check list

(see reverse for more details)

 Extended Warranty must be purchased prior to the expiration of active warranty







Maintenance Checklist

Our comprehensive 26 point maintenance checklist are thorough, ensuring the Automet is running efficiently.

26 Point Maintenance Checklist Power Head

- Theck head rotation for smooth operation
- $oxed{oxed}$ Check and adjust belt tension
- ☐ Check proper pressure of single & central load
- ☑ Check head post rotation and lubricate if necessary.
- Check upper and lower limit switches for proper head setting
- ☑ Check and verify platen function properly
- ✓ Replace bearings
- ✓ Monitor motor performance
- ☑ Replace platen screws and o-rings
- ☑ Check cooling fan and clean fan filter
- $oxedsymbol{oxed}$ Check and verify operation of machine.
- ✓ Check for proper operation of the waterspout retracting tube.

- ☑ Check head movement for smooth operation
- M Check all air lines for leaks
- ☑ Check head LED lights
- ☑ Ensure proper function of head controls (start buttons)

Base

- ☑ Check and clean water drain hole for proper drainage
- ☑ Replace solenoid, water filter, and water valve
- ✓ Check all water lines for leaks.
- ☑ Replace bowl liner and clean bowl
- ☑ Replace and adjust belt tension
- ☑ Update firmware to latest release



Wear Parts Kit

The parts listed below are included with every Preventative Maintenance visit


AutoMet Wear Parts Kit 49725SRV

Includes

Platen O-Rings Water Solenoid Water Valve Water Filter Kit

Platen Bolt

Disposable Bowl Liners Platen Drive Belt 250 main bearings Platen Bearings Kit Intake Water Hose

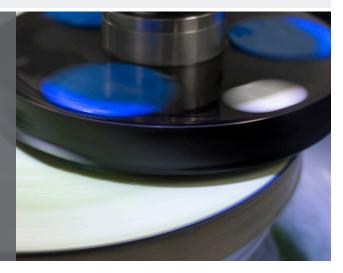



BUEHLER Worldwide Headquarters North America-South America Offices

41 Waukegan Road Lake Bluff, Illinois 60044-1699 USA P: 800 BUEHLER (800-283-4537) P: (847) 295-6500 www.buehler.com | info@buehler.com European Headquarters BUEHLER Germany info.eu@buehler.com

BUEHLER France info.eu@buehler.com

BUEHLER United Kingdom info.eu@buehler.com

BUEHLER China info.cn@buehler.com

BUEHLER Japan info.japan@buehler.com

BUEHLER Asia-Pacific info.asia@buehler.com